

Protocol	SCHOTT	
NEXTERION® IC-16 and IC-16s Cleaning protocol	Dok-Nr.:	LS6-HBM-M-002
	Version:	1.0
	Seite:	1/3
	Datum:	© June 2011

1	Cleaning after usage with non-biohazardous components:.....	2
2	Cleaning after usage with biohazardous components:.....	3

For Technical Assistance, please contact

**SCHOTT Technical Glass
Solutions GmbH**

Otto-Schott-Straße 13
07745 Jena
Germany

Phone: +49-(0)3641-681-4069

Fax: +49-(0)3641-681-4970


E-Mail: coatedsubstrate@schott.com E-Mail:
coatedsubstrate@schott.com

Additional information at:
www.schott.com/nexterion

Protocol	SCHOTT	
NEXTERION® IC-16 and IC-16s Cleaning protocol	Dok-Nr.:	LS6-HBM-M-002
	Version:	1.0
	Seite:	2/3
	Datum:	© June 2011

1 Cleaning after usage with non-biohazardous components

- 1.1 After usage immediately rinse the disassembled IC-16 chamber and superstructure with deionised water.
- 1.2 Submerge all components in a vessel containing Liquinox™ or a similar cleaning solution.
- 1.3 Soak for 30 minutes or longer.
- 1.4 Using a nylon brush scrub all components taking care to ensure the bristles scrub the inside of the wells on the superstructures and upper plate.
- 1.5 Rinse with extensive amounts of deionised water.
- 1.6 Rinse with extensive amounts of double-distilled/Millipore™ water.

Protocol		
NEXTERION® IC-16 and IC-16s Cleaning protocol	Dok-Nr.:	LS6-HBM-M-002
	Version:	1.0
	Seite:	3/3
	Datum:	© June 2011

2 Cleaning after usage with biohazardous components

- 1.1 After usage immediately submerge the disassembled IC-16 chamber and superstructure in a vessel containing a $\geq 10\%$ bleach solution ($\geq \sim 0.5\%$ sodium hypochlorite).
- 1.2 Soak for 30 minutes or longer.
- 1.3 Rinse all components with extensive amounts of deionised water.
- 1.4 Submerge the superstructures in a vessel containing 70 % Ethanol.
- 1.5 Soak for a 1 hour to 24 hours.
- 1.6 Rinse the superstructures in extensive amounts of deionised water.
- 1.7 Submerge all components in a vessel containing Liquinox™ or a similar cleaning solution.
- 1.8 Soak for 30 minutes.
- 1.9 Using a nylon brush scrub all components taking care to ensure the bristles scrub the inside of the wells on the superstructures and upper plate.
- 1.10 Rinse with extensive amounts of deionised water.
- 1.11 Rinse with extensive amounts of double-distilled/Millipore™ water.
- 1.12 Tap the chamber pieces gently to remove excess liquid.
- 1.13 Place chamber and superstructure on a paper towel to dry.
- 1.14 Dispose any superstructures which have abrasions or roughened surfaces.